

## THE ENVIRONMENTAL COMPLIANCE EXPERTS



| DESCRIPTION  | QTY |
|--|-----|
| FACE MASK & EYE PROTECTION   | 5   |
| PROTECTIVE APRON   | 5   |
| 500G PREMIUM BODY FLUID ABSORBENT (INCLUDING ZEOLITE WITH ADDED ODOUR CONTROL) | 2   |
| SCOOP & SCRAPER  | 5   |
| YELLOW CLINICAL WASTE BAG  | 10  |
| CABLE TIE  | 10  |
| PH NEUTRAL DETERGENT   | 2   |
| ABSORBENT CLEANING CLOTH   | 10  |
| ANTISEPTIC TOWELETTE   | 10  |

## BIOHAZARD SPILL CLEANUP PROCEDURE

- 1 Prepare:** Open the Biohazard Spill Kit and lay out contents on a clean surface.
- 2 Wear PPE:** Put on the Face Mask, Eye Protection, Apron, and Gloves.
- 3 Apply Absorbent:** Evenly pour BioSmart Zeolite Powder over the spill, working from the edges inward.
- 4 Absorb the Spill:** Ensure full coverage and stir with the scraper if needed until liquid is absorbed.
- 5 Dispose of Waste:** Use the Scoop & Scraper to transfer waste into a yellow Clinical Waste Bag. Seal with a cable tie and double-bag using the second yellow bag.
- 6 Clean Residue:** Use BioSmart pH Neutral Detergent and a cloth to wipe the area. Dispose of the cloths in the second yellow waste bag.
- 7 Remove PPE:** Place all used PPE (mask, gloves, apron, eye protection) into the second yellow waste bag and seal with a cable tie.
- 8 Sanitize Hands:** Wash with soap and water or use antiseptic wipes if no sink is available.
- 9 Dispose Properly:** Follow Local Government Regulations for biohazard waste disposal.

**To re-order call 0434 314 803**